Safety Instructions PM2 Photometer

PENDOTECH
Leading Process Analytics

Contents

en	Safety Instructions	3
de	Sicherheitshinweise	6
fr	Consignes de sécurité	9
es	Instrucciones de seguridad	12
it	Istruzioni sulla sicurezza	15
zh	安全说明	18
cs	Pokyny k bezpečnosti	20
da	Sikkerhedsvejledning	23
et	Ohutustehnika juhend	26
fi	Turvaohjeet	29
el	Οδηγίες ασφαλείας	32
hu	Biztonsági előírások	35
ga	Treoracha um Shábháilteacht	38
ja	安全ガイド	41
ko	안전 지침	44
lt	Drošības instrukcijas	47
lv	Saugaus darbo instrukcijas	50
mt	Struzzjonijiet tas-Sigurtà	53
nl	Veiligheidsvoorschriften	56
pl	Instrukcje dotyczące bezpieczeństwa	59
pt	Instruções de Segurança	62
ru	Правила техники безопасности	65
sv	Säkerhetsinstruktioner	68
sl	Varnostna navodila	71
sk	Pokyny k bezpečnosti	74

METTLER TOLEDO

en Safety Instructions

1 Definition of Equipment and Documentation Symbols and Designations

WARNING: POTENTIAL FOR PERSONAL INJURY

CAUTION: Possible instrument damage or malfunction

NOTE: Important operating information

On the photometer or in the manual text indicates: Caution and/or other possible hazard including risk of electric shock (refer to accompanying documents).

2 Each prospective user must test the measurement unit for its proposed application to determine its suitability for the purpose intended prior to incorporating the sensor to any process or application. Proper safeguards must be put into place for the process in which the unit is used.

The functioning and operational safety of the product can only be ensured if the user observes the usual safety precautions as well as the specific safety guidelines stated in these operating guidelines:

3 Warnings

- The PM2 Photometer must only be operated under the specified operating conditions.
- The PM2 Photometer should be installed and operated only by personnel familiar with the Photometer and who are qualified for such work.
- Before connecting the device to a supply unit, ensure its output voltage cannot exceed 48 VDC. Do not use alternating current or a main power supply!
- **Warning:** Only use a power supply with double insulation or reinforced insulation to supply this product.
- **Warning:** Do not submerge this product. Protect the product before cleaning with any liquids by covering openings that expose the internal components.
- Except for routine maintenance and cleaning procedures the PM2 Photometer must not be tampered with or altered in any manner.

3 Warnings

- Maintenance, and repair work must only be carried out by PendoTECH.
- METTLER TOLEDO/PendoTECH accepts no responsibility for damage caused by unauthorized modifications to the Photometer.
- Follow all warnings, cautions, and instructions indicated on and supplied with this product.
- Install equipment as specified in this instruction manual. Follow appropriate local and national codes.
- Electrical installation must be per the National Electrical Code and/or any other applicable national or local codes.
- Power supply must be disconnected before servicing.
- If this equipment is used in a manner not specified by the manufacturer, its protection against hazards may be impaired.
- The safety of the user may be endangered if the instrument:
 - is visibly damaged
 - no longer operates as specified
 - has been damaged in transport
- Relay action: the PM2 Photometer relays will always de-energize on the loss of power, equivalent to a normal state, regardless of the relay state setting for powered operation. Configure any control system using these relays with fail-safe logic accordingly.
- Process upsets: because process and safety conditions may depend on consistent operation of this photometer, provide appropriate means to maintain operation during sensor cleaning, replacement or sensor or instrument calibration.
- The terminals are suitable for single wires / flexible leads 0.2 mm² to 1.5 mm² (AWG 16 – 24).

4 **Optical Safety**

- **Warning:** The photometer can emit light in the range of 200 nm – 1100 nm. Hazardous UV and IR radiation are emitted from the unit. Never directly stare into the source/return ports on the photometer or the fiber optic cables, as this can cause eye damage. The viewer-related risk depends on how the users install and use the product.
- Ensure that both ends of the fiber optic cables (photometer and flow cell connections) are properly secured (firmly hand-tight) before powering on the photometer.

-
- The photometer should always be powered off when handling the fiber optic cables. Protective eyewear must be worn if you must handle the fibers when the unit is powered on.
 - In the event of product failure, do not attempt to open the unit or replace the LED. There are no user-serviceable parts.

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All rights reserved